

OBTAIN A MORE COMPLETE PROFILE OF SAMPLE COMPONENTS AND THE FREEDOM TO USE UV-INTERFERING MOBILE PHASES

Application note # QAN-0073

PARAMETERS

NQAD


GAS: Air
FILTER: 1.3 sec
GAIN: 1X
EVAPORATION TEMP: OFF

HPLC

MOBILE PHASE A: 5 mM ammonium acetate
MOBILE PHASE B: 0.05% formic acid in ACN
METHOD: 25:75 A:B

FLOW RATE: 0.5 mL/min
COLUMN: Tosoh amide 80, 2.0 x 150 mm, 5u
COLUMN TEMP: Ambient
INJECTION VOLUME : 10 µL
ANALYTE: Melamine
CONCENTRATION: 100 ng

CHROMATOGRAMS

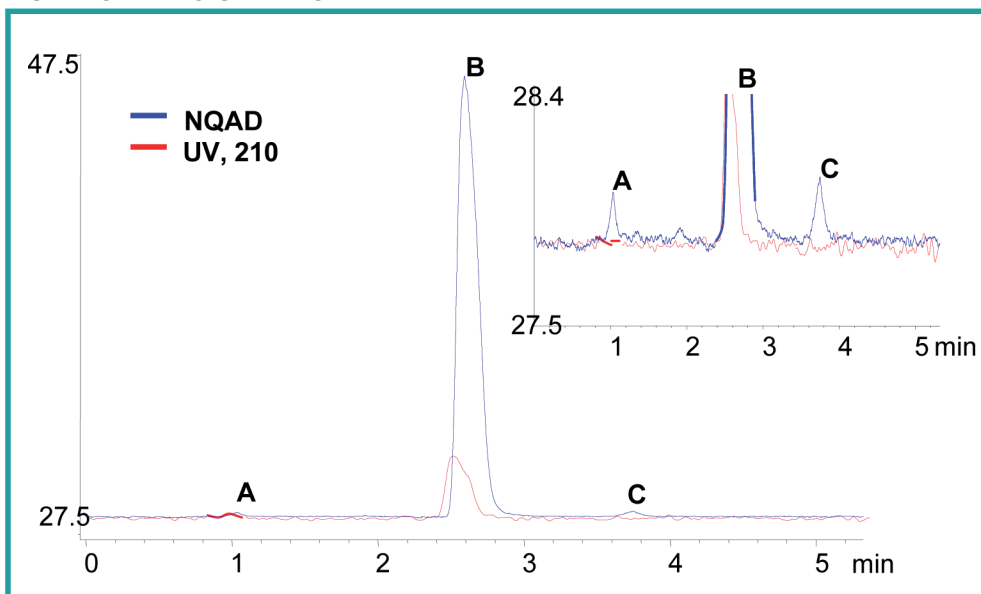

(A) = cyanuric acid, (B) = melamine, (C) = unidentified impurity



ERC GmbH
Otto-Hahn-Strasse 28-30
D-85521 Riemerling
☎ +49(0)89/660 55 696
☎ +49(0)89/608 24 826
partner@erc-hplc.de
www.erc-hplc.de